



VERSATILE, HIGH GRADE DISINFECTION WITH EASY APPLICATION



Klercide™ 70|30 Pharma Ethanol blended with WFI

70|30 Pharma Ethanol denatured with Isopropyl Alcohol blended with Water for Injection (WFI)

The solution of ethanol with isopropyl alcohol as the denaturant will help ensure that residues on evaporation are comparable to USP grade ethanol.

Effective for use in direct product contact areas, the product is formulated without non-volatile denaturants, leaving no surface residue contaminants.

Klercide 70|30 Pharma Ethanol is:

- ▲ 0.2 micron filtered, filled and bagged in a Grade B cleanroom
- ▲ Available double and triple bagged
- ▲ Equipped with the SteriShield Delivery System

The SteriShield Delivery System (SDS) features:

- ▲ A closed trigger system providing a completely sealed system
- ▲ Validated in-use shelf life of three months
- ▲ A fully adjustable spray pattern to ensure optimum surface coverage

PRODUCT QUALITY ASSURANCE

Manufactured in our state-of-the-art facilities to the principles of cGMP, Klercide 70|30 Pharma Ethanol is safe for use in critical aseptic core manufacturing areas.



Klercide™ 70|30 Pharma Ethanol

Blended with WFI



70|30 Pharma Ethanol denatured with Isopropyl Alcohol blended with Water for Injection (WFI)

Why use Klercide 70|30 Pharma Ethanol?

EFFICIENT

Minimal residues and optimum surface coverage ensure the spray covers large areas effectively, while not leaving residues that may create risk in critical areas. The trigger spray can be adjusted to provide a jet or a fine spray to ensure thorough wetting of the surface, so that disinfection procedures are effective.

COST EFFECTIVE

Klercide 70|30 Pharma Ethanol can be completely dispensed from the bottle so there is no wastage. The trigger spray uses the SteriShield Delivery System which has been validated to enable an in-use shelf life of three months, so there is no need to discard unused product, therefore reducing waste.

DESIGNED FOR CLEANROOMS

Klercide 70|30 Pharma Ethanol is manufactured in Ecolab's ISO 9001 accredited facilities to the principals of cGMP. These manufacturing conditions ensure Klercide 70|30 Pharma Ethanol is manufactured to the highest quality and are compliant for use in critical areas.

Klercide 70|30 Pharma Ethanol is:

- ▲ Endotoxin tested to an approved pharmacopoeia method
- ▲ Supported with Certificates of Analysis and Sterilization with every batch
- ▲ Supported with full documentation

MICROBIOLOGICAL TEST RESULTS

EFFICACY	CONTACT TIME
Disinfection	5 min
Virucidal	5 min
Fungicidal	10 min and 5 min*
Non-Food Contact Sanitizer	3 min

*Please see label for specific claim

WORLDWIDE HEADQUARTERS

1 Ecolab Place
St. Paul, MN 55102
USA

www.ecolab.com/lifesciences
www.ecolab.com/pharmaceutical

© 2019 Ecolab USA Inc. All rights reserved. 30JAN19/NA

PHARMACEUTICAL

Efficient. Simple. Economical.

APPLICATION AREA

For use on floors, walls and hard surfaces.

INSTRUCTIONS FOR USE

Spray: For optimum results hold bottle upright and spray from a distance of 6-8 inches. Always close the nozzle after use.

Wetting: For optimum results ensure complete wetting of the surfaces using the spray and wipe technique.

INGREDIENTS

70% ethanol denatured with isopropyl alcohol and blended with Water for Injection.

REGISTRATION DETAILS

EPA Reg. No. 1677-252
Canadian DIN 02453002

PACKAGING INFORMATION

Klercide 70|30 Pharma Ethanol

UNIT SIZE	UNIT OF SALE
32 oz spray	6 x 32 oz.

FOR PRODUCT ORDER CODES AND ADDITIONAL INFORMATION, PLEASE CONTACT YOUR ACCOUNT MANAGER OR VISIT OUR WEBSITE

SHELF LIFE

Unopened, this product has a three year shelf life.

ALWAYS READ THE LABEL AND PRODUCT INFORMATION BEFORE USE.

ECOLAB®